

FlexAble CoraLite® Plus 405 Antibody Labeling Kit for Rabbit IgG

Catalog Number: KFA006

Product Information

| | |
|--|--|
| Product name | FlexAble CoraLite® Plus 405 Antibody Labeling Kit for Rabbit IgG |
| Assay type | Antibody labeling |
| Tested applications | IF, FC |
| Species Reactivity | Rabbit IgG |
| Antibody amount per labeling reaction | 0.5 µg antibody |
| Conjugate | CoraLite® Plus 405 |
| Excitation / Emission maxima wavelengths | 399 nm / 422 nm |

Kit Components

| Component | 10 rxns | 50 rxns | 4×50 rxns |
|--|---------|---------|-----------|
| CoraLite® Plus 405 FlexLinker for Rabbit IgG | 10 µL | 50 µL | 4×50 µL |
| FlexQuencher for Rabbit IgG | 20 µL | 100 µL | 4×100 µL |
| FlexBuffer | 100 µL | 500 µL | 4×500 µL |

包装规格

10/50/4×50 reactions

Storage Condition

Store for 1 year at -20°C or for 6 months at +4°C upon receipt. Avoid exposure to light.

FAQ

Q: What are the FlexLinker, FlexQuencher and FlexBuffer?

A: The FlexLinker is a small polypeptide to which dyes are covalently conjugated that can label unconjugated primary antibodies. The FlexQuencher is an Fc-containing fragment that neutralizes the excess FlexLinker. The FlexBuffer is a PBS-based buffer.

Q: What is the largest quantity I can label?

A: With a standard kit size (50 reactions), you can label 25 µg of one antibody or up to 50 different antibodies. You can easily scale up the antibody amount per labeling approach.

Q: What is the lowest concentration of my primary antibody that I can use?

A: Our protocol uses 0.5 µg of primary antibody in 7 µL, which ends up at 0.07 mg/mL. If the concentration of your antibody is lower, you can also use a larger volume than 7 µL.

Q: Can I label primary antibodies stored in BSA, glycerol, Tris buffer and/or preservatives?

A: Yes, FlexAble Antibody Labeling Kits have been validated with carriers and amine buffers. Neither BSA nor amine buffers, in any chosen concentration, interfere with the labeling. 50% glycerol as well as preservatives like sodium azide are also compatible with the kit.

Q: How many different primary antibodies can I label with one kit?


A: You can label up to 50 different antibodies with our FlexAble 50 rxn Kit, and up to 10 antibodies with our FlexAble 10 rxn Kit.

Q: Will I observe cross-reactivity/leaking when I use two FlexAble-labeled antibodies from the same species during multiplexing?


A: FlexAble labels primary antibodies with a high affinity FlexLinker. Dissociation of FlexLinker from one antibody and association to another antibody is rare. If you observe leaking, we recommend adding more FlexQuencher to remove unbound FlexLinker, or you can try sequential staining of the labeled antibodies.

[More FAQs](#)


Validation Data

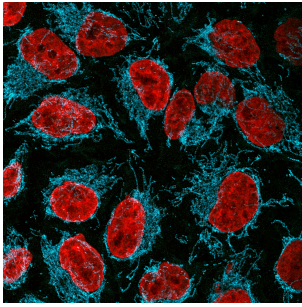

Immunofluorescence of HeLa: PFA-fixed HeLa cells were stained with anti-TOM70 (14528-1-AP) labeled with FlexAble CoraLite® Plus 405 Kit (KFA006, cyan). Cell nuclei are in red. Confocal images were acquired with a 100x oil objective and post-processed. Images were recorded at the Core Facility Bioimaging at the Biomedical Center, LMU... Munich.

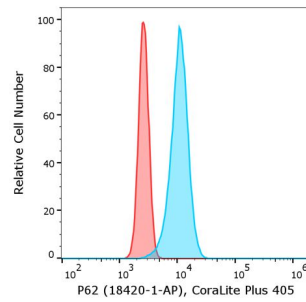

Flow cytometry of PBMC. 1×10^6 human peripheral blood mononuclear cells (PBMCs) were intracellularly stained with 0.5 μ g anti-human P62 (SQSTM1) antibody (18420-1-AP) labeled with FlexAble CoraLite® Plus 405 Kit (KFA006, cyan) or with 0.5 μ g isotype control antibody labeled with FlexAble CoraLite® Plus 405 Kit (KFA006, red).

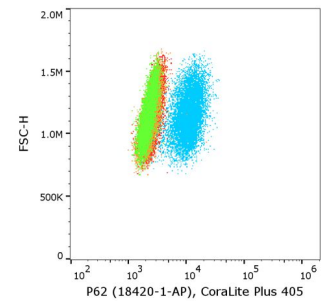

Flow cytometry of HEK293. 1×10^6 HEK293 cells were intracellularly stained with 0.5 μ g anti-human P62 (SQSTM1) antibody (18420-1-AP) labeled with FlexAble CoraLite® Plus 405 Kit (KFA006, blue) or with 0.5 μ g isotype control antibody labeled with FlexAble CoraLite® Plus 405 Kit (KFA006, red) or with FlexLinker CoraLite® Plus 405 for rabbit... IgG (orange). Unstained HEK293 cells were used as a reference (red).

For technical support and original validation data for this product please contact

T: 027-87531629

E: Proteintech-CN@ptglab.com

W: ptgcn.com

This product is exclusively available under Proteintech Group brand and is not available to purchase from any other manufacturer.